

CULTURING OF *PSEUDO-NITZSCHIA* SPECIES ON AGAR-F/2 MEDIUM

PREAMBLE. The method outlined below was employed for *Pseudo-nitzschia* species. The long-term cultivation of this pennate diatom in liquid media is difficult and often unsuccessful. The growth on agar plates can be considered an alternative method which ensures a reduction of operational time and good results in terms of survival/shape of the cultures.

1. Preparing of 1 l of agar-f/2 solution (at 0.2% of agar)

Add 2 gr of agar to 1 l of seawater.

Sterilize by autoclave.

Add stock solutions as reported in the recipe for F/2 medium in Guillard (1975).

N.B. Add the stock solutions when the agar has still not reached the gelling point.

Prepare 6-wells plates with 4 ml of agar-f/2 at 0.2%.

Bring the plates at room temperature, seal them with a strip of Parafilm™ to minimise dehydration of the agar. The cultures should be inoculated the next day.

2. Plating

a) From cultures maintained in liquid medium to agar:

Select a dense, healthy culture in exponential growth phase. Add 0.2 ml of the culture and 1ml of f/2 medium to an agar-well and mix it with a loop to be sure that the cells are homogeneously mixed into the agar.

b) From agar to agar:

Add 0.5 ml of f/2 medium to the culture on agar to be transferred and mix it with a loop to homogenise it. In this way, the agar medium becomes softer and favours the success of culture transfer.

Isabella Percopo, Carmen Minucci,
Marina Montresor, Diana Sarno
Stazione Zoologica Anton Dohrn
Villa Comunale 80121 - Napoli (Italy)

*Algae: Method has been successfully
employed for Pseudo-nitzschia species*

Apparatus:

Solutions: F/2 medium, agar

Plasticware: 6-wells plates, loops

*Additional information: Guillard, R.R.L.
1975. Culture of phytoplankton for
feeding marine invertebrates. pp 26-
60. In Smith W.L. and Chanley M.H
(Eds.) Culture of Marine Invertebrate
Animals. Plenum Press, New York, USA.*

ASSEMBLE
ASSOCIATION OF EUROPEAN MARINE BIOLOGICAL LABORATORIES

Using the loop pick some material (microalgae and agar) from the old plate and mix it in the new agar-well.

Repeat this step for 4-5 times for the same culture strain, to be sure that enough material is transferred in the agar-well.

Add 1 ml of f/2 medium and mix again.

Seal the plate with a strip of Parafilm™.

Keep the cultures at 12:12 L:D cycle; 18-20°C and 50 $\mu\text{mol photons m}^{-2} \text{sec}^{-1}$ of irradiance. The temperature can vary depending from the sites where strains have been isolated.